

MONOTOPE™ product information for

Plasmodium falciparum

I. Monoclonal Antibody (Mouse) Specificity

Product No.'s	Ig class	
3564 NA	IgG2a	Recognizes histidine rich protein II of P. falciparum. ELISA. See pairing recommendations below.
3567 NA	IgG2a	Recognizes histidine rich protein II of P. falciparum. ELISA. See pairing recommendations below.
3568 NA	IgG2b	Recognizes histidine rich protein II of P. falciparum. ELISA. See pairing recommendations below.
3569 NA	IgG2b	Recognizes histidine rich protein II of P. falciparum. ELISA. See pairing recommendations below.
3566 NA	IgG1	Recognizes merozoite surface protein 1 (MSP1) of P. falciparum. ELISA & WB.

II. Purified Preparations

Product No.'s	3564
	3567
	3568
	3569
	3566

MONOTOPE™ purified preparations consist of >90% pure mouse monoclonal antibody which has been purified from ascites fluid or culture medium by protein A chromatography or sequential differential precipitations. The final preparation is formulated to a protein concentration of 100 µg/ml in 0.01 M phosphate buffered saline, pH 7.2 and contains 0.1% sodium azide. Each vial contains 1.0 ml. This product contains no stabilizing proteins and should be stored at 2-8°C until ready for use. Working dilution must be determined by the user. Suggested starting ranges are 1:10-1:50 for IFA and 1:20-1:200 for ELISA.

THESE PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES

III. Fluorescein Conjugates

Product No.'s	NA
	NA
	NA
	NA
	NA

These MONOTOPE™ products consist of purified monoclonal antibody conjugated with high purity isomer I of fluorescein isothiocyanate. Care is taken to ensure complete removal of any free fluorescein from the final product. The final preparation is formulated to an antibody concentration of 100 µg/ml in 0.01 M phosphate buffered saline, pH 7.2 containing 0.1% sodium azide plus bovine serum albumin at 10 mg/ml. Each vial contains 1.0 ml. This product should be stored at 2-8°C until ready for use. Avoid repeated freeze-thawing by storing multiple aliquots at -20°C. Applications for these products include direct FA staining of target antigen in a permissive tissue culture system. Working strength must be determined by the user for each specific application but a starting range of 1:5 - 1:20 is recommended. Acetone fixation of the antigen source is recommended prior to staining.

THESE PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES

Comments: Best ELISA pairs (capture - conjugate):

3564 - 3567
3568 - 3564
3569 - 3564