

MONOTOPE™ product information for

Mycoplasma pneumoniae P1

I. Monoclonal Antibody (Mouse) Specificity

Product No.'s	Ig class	
6532 NA	IgG2b	Specific for the P1 adhesin of M. pneumoniae. ELISA & IFA. See pairing recommendations below.
6533 NA	IgG1	Specific for the P1 adhesin of M. pneumoniae. ELISA & IFA. See pairing recommendations below.
6534 NA	IgG2b	Specific for the P1 adhesin of M. pneumoniae. ELISA & IFA. See pairing recommendations below.
6535 NA	IgG2b	Specific for the P1 adhesin of M. pneumoniae. ELISA & IFA. See pairing recommendations below.
6536 NA	IgG2a	Specific for the P1 adhesin of M. pneumoniae. ELISA & IFA. See pairing recommendations below.
6537 NA	IgG2a	Specific for the P1 adhesin of M. pneumoniae. ELISA & IFA. See pairing recommendations below.

II. Purified Preparations

Product No.'s	6532
	6533
	6534
	6535
	6536
	6537

MONOTOPE™ purified preparations consist of >90% pure mouse monoclonal antibody which has been purified from ascites fluid or culture medium by protein A chromatography or sequential differential precipitations. The final preparation is formulated to a protein concentration of 100 µg/ml in 0.01 M phosphate buffered saline, pH 7.2 and contains 0.1% sodium azide. Each vial contains 1.0 ml. This product contains no stabilizing proteins and should be stored at 2-8°C until ready for use. Working dilution must be determined by the user. Suggested starting ranges are 1:10-1:50 for IFA and 1:20-1:200 for ELISA.

THESE PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES

III. Fluorescein Conjugates

Product No.'s	NA
	NA
	NA
	NA
	NA
	NA

These MONOTOPE™ products consist of purified monoclonal antibody conjugated with high purity isomer I of fluorescein isothiocyanate. Care is taken to ensure complete removal of any free fluorescein from the final product. The final preparation is formulated to an antibody concentration of 100 µg/ml in 0.01 M phosphate buffered saline, pH 7.2 containing 0.1% sodium azide plus bovine serum albumin at 10 mg/ml. Each vial contains 1.0 ml. This product should be stored at 2-8°C until ready for use. Avoid repeated freeze-thawing by storing multiple aliquots at -20°C. Applications for these products include direct FA staining of target antigen in a permissive tissue culture system. Working strength must be determined by the user for each specific application but a starting range of 1:5 - 1:20 is recommended. Acetone fixation of the antigen source is recommended prior to staining.

THESE PRODUCTS ARE FOR RESEARCH USE ONLY. NOT FOR USE IN DIAGNOSTIC PROCEDURES

Comments: Best ELISA pairs (capture - conjugate): 6532-6535, 6533-6535, 6534-6537, 6533-6536

These antibodies do not cross react with: Ureaplasma urealyticum, Mycoplasma hominis or Mycoplasma genitalium